


interpretation

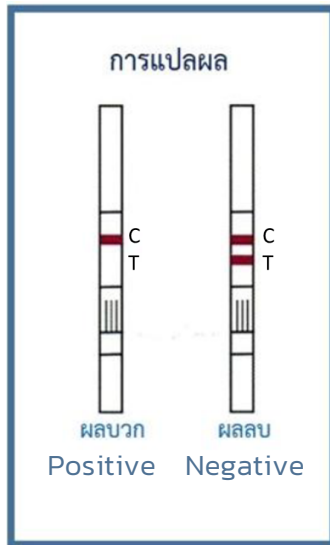

- **Positive:** 1 line (appear at the "C" area) means THC concentration is higher than 0.2%.
- **Negative:** 2 lines (appear at the "C" and "T" areas) mean THC concentration is lower than 0.2%.

Note: If the line at "C" does not appear, the result is invalid.



Department of Medical Sciences
Ministry of Public Health, Thailand

Contact us

88/7 Tiwanon Rd. Amphoe Muang,
Nonthaburi 11000
www.dmsc.moph.go.th



Test Kann

Test kit for THC in
cannabis extract and oil



Test kit for THC
in cannabis extract and oil



THC test kit

Notification of the Ministry of Public Health Re: Specification of Type 5 Narcotics B. E. 2565 (2022) has prescribed that "the extract from every part of cannabis and hemp which contains tetrahydrocannabinol (THC) higher than 0.2% is specified as type 5 Narcotics."

Test samples


"Cannabis extract and oil (to test whether it contains THC > 0.2%)



Test kit components



1. Testing cassette
2. solution 1
3. solution 2
4. plastic spoon
5. dropper

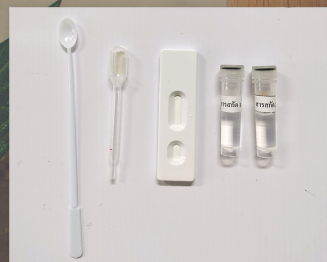

Test procedure




1




Measure 1 spoon of cannabis extract or oil by using the plastic spoon.

2

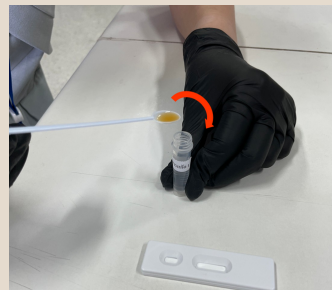


Pour the sample from the spoon into the tube of solution 1.

3

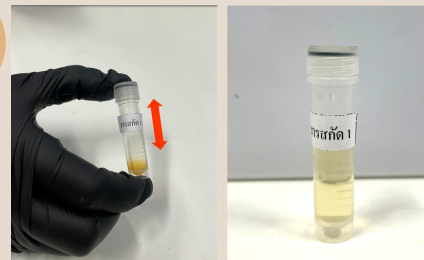


Vertically shake the tube to mix it thoroughly.

4

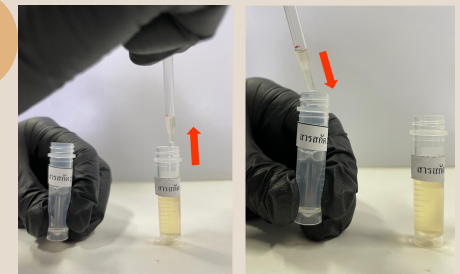


Use the dropper to suck solution 1 up to the specified level and drop it into the solution 2 tube.

5

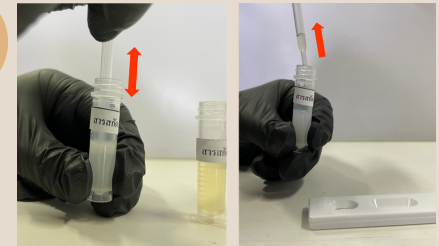

Mix solution 2 in the tube by using dropper then suck the solution up to the specified level.

6

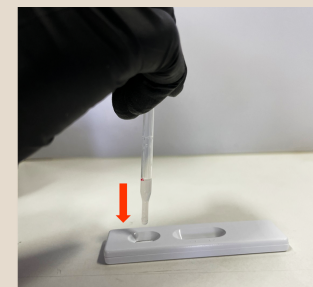

Drop the solution into the sample well of the testing cassette, wait for 15 min, and read the result.